



# How to Install An USB Key

1. Double click the exe data "WkRuntime.exe" and follow the direction on the screen.  
 ※ Do not insert the USB key to your PC yet at this moment.


① Double click "WkRuntime.exe"

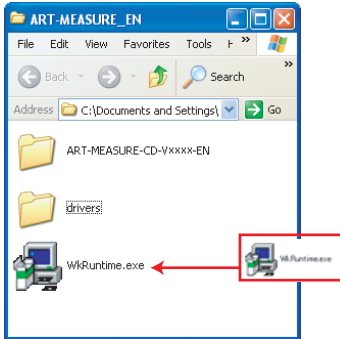


② Click "NEXT"

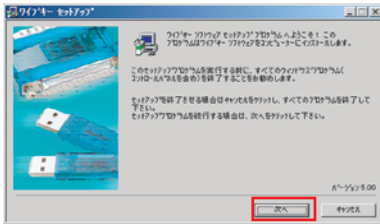


③ Choose a language

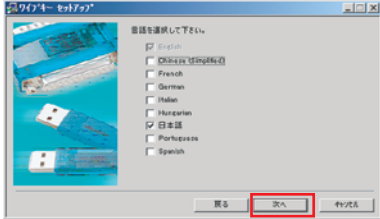


④ Click "NEXT"

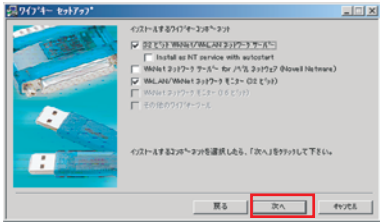



⑤ Click "NEXT"

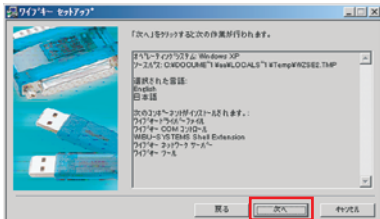

⑥ Click "NEXT"

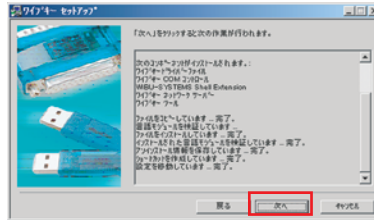


⑦ Click "FINISH"

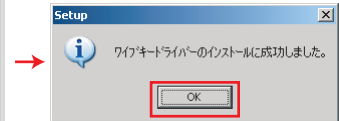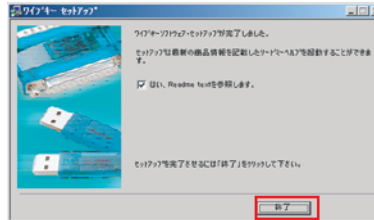

2. Insert the USB key to your PC.



⑧ Choose 「Install the software automatically [Recommended]」. And click "NEXT".

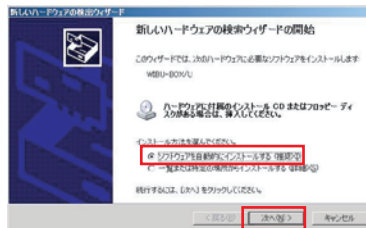

USB Key

The screen on left is shown.

⑨ Start installing.

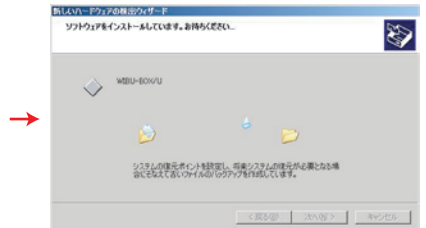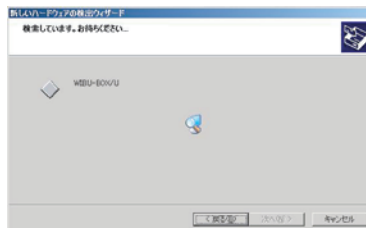

⑩ Click "FINISH" to finish installation.



3. After install the measure software (See next page for more detail), start the software.  
 \*You cannot use the software without USB key.